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## Instruments and Software

Item	Description	Part No.	Unit
Oven	ChroZen GC Mainframe Assembly with UPC Detector Board Unit	6701012502	1
Inlet	Capillary Inlet Assembly for ChroZen GC (毛细管)	6701012550	1
	Packed Inlet Assembly for ChroZen GC (填充柱)	6701012510	2
Detector	FID Assembly for ChroZen GC	6701012590	1
	TCD (Thermal Conductivity Detector) Assembly for ChroZen GC	6701012570	2
CDS	YL-Clarity software for single instrument of YCM GC	5301011020	1
Column	GS-gaspro 60 m 0.32 mm	-	1
	2m 1/8 2mm MOLECULAR SIEVE 5A 60/80	-	2
	6Ft Porapak Q 60/80 SS	-	2
Install. Option	Start-up kit	1601011110	1
Valve	Automatic Gas Valve, 2 pos/6 port, Valcon E/Max: 225°C for ChroZen GC Micro-electric actuator type with 250 ul sample loop	6501011250	2
	Automatic Gas Valve, 2 pos/10 port, Valcon E/Max: 225°C for ChroZen GC Micro-electric actuator type with 250 ul sample loop	6501011280	2

Fig 1. ChroZen GC with Column Switching Valve System

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Initial



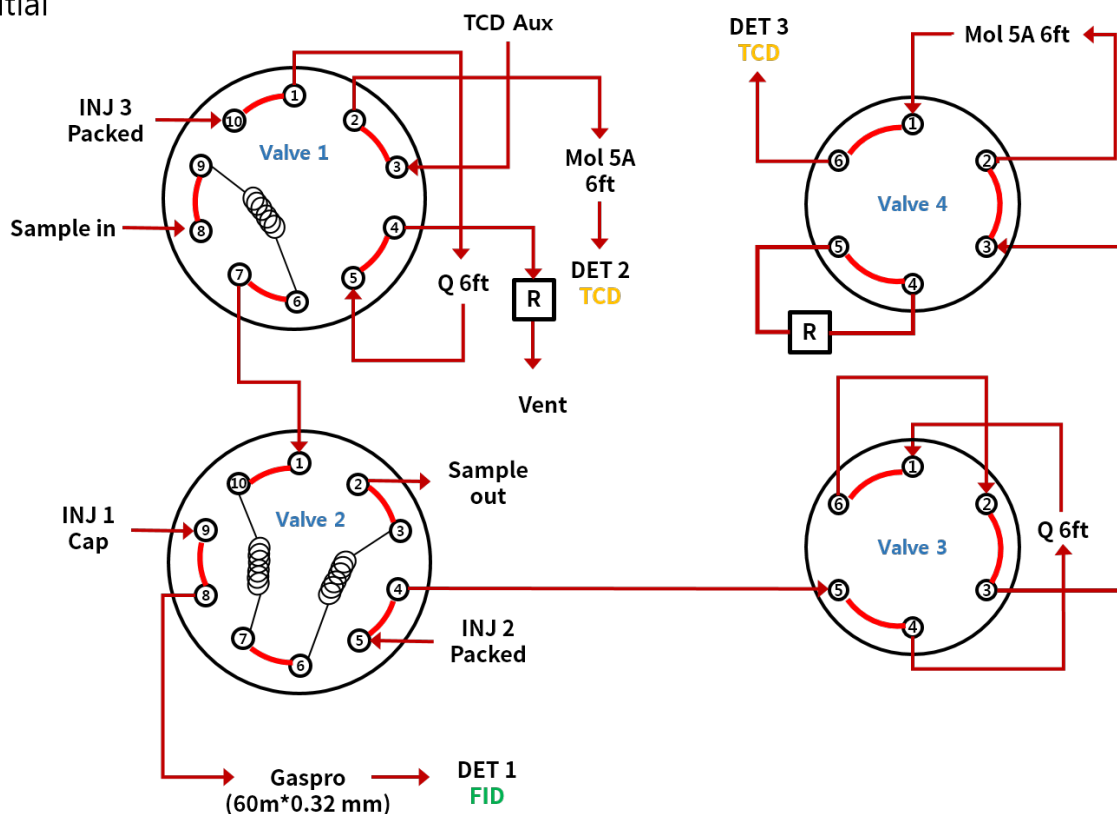
Fig 2. Valve Diagram of Natural Gas Analysis

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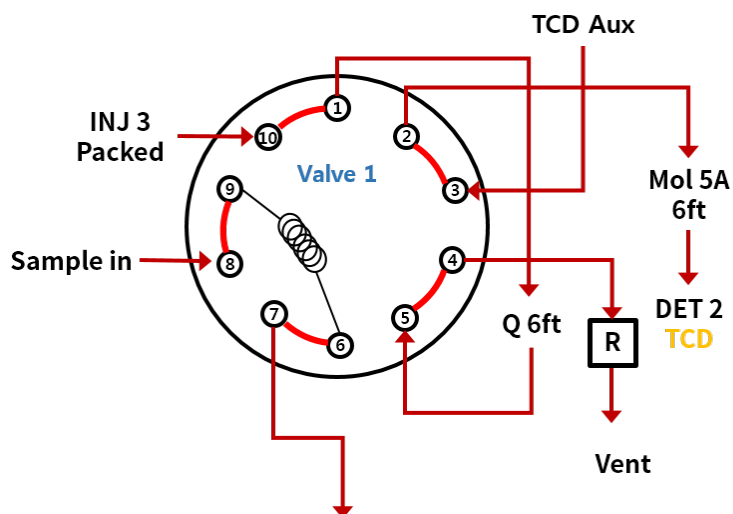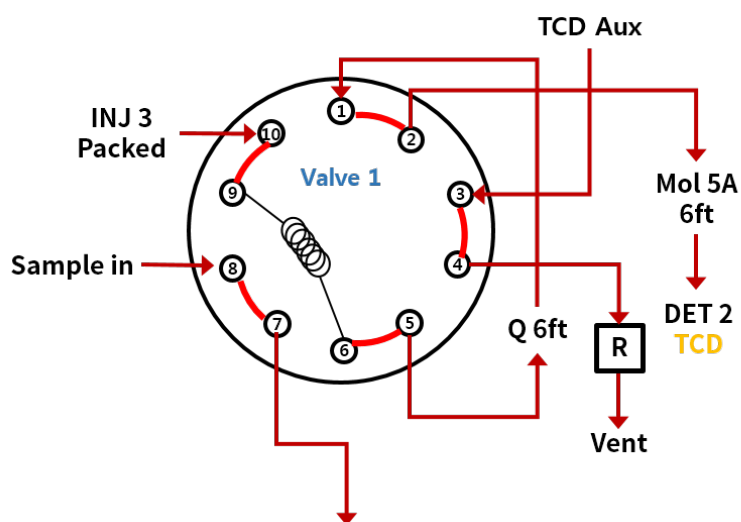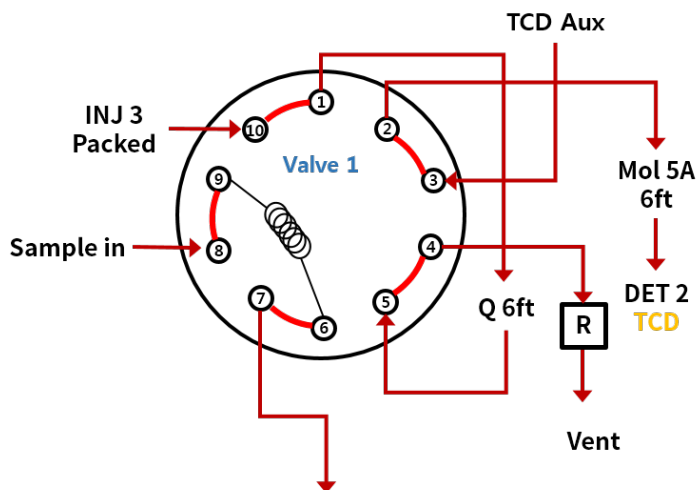
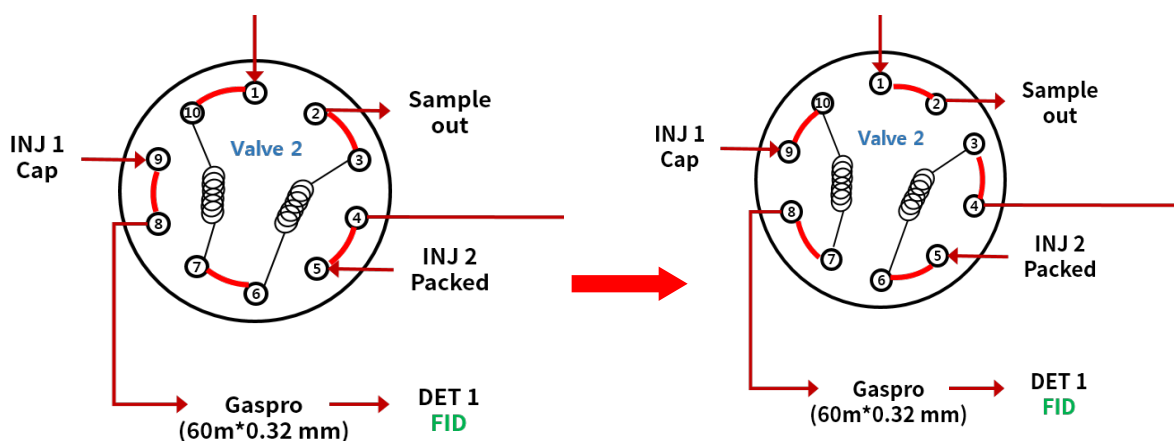
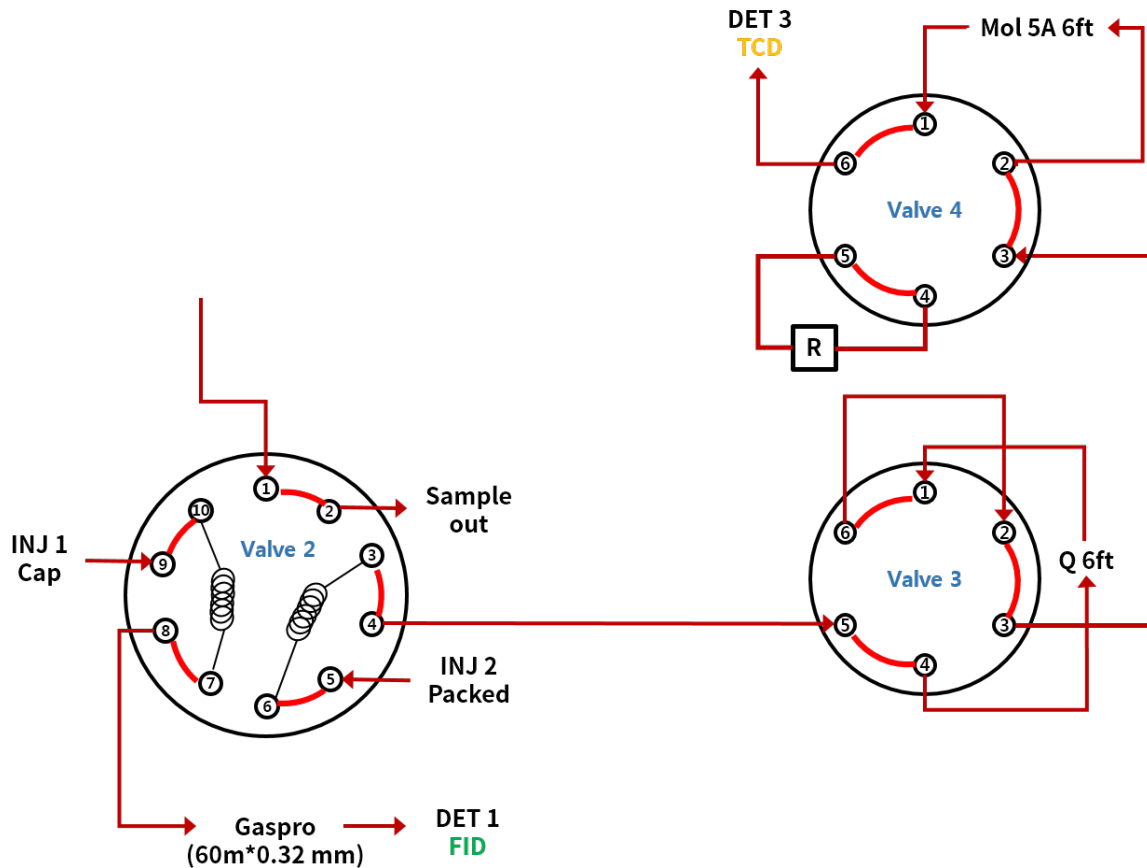
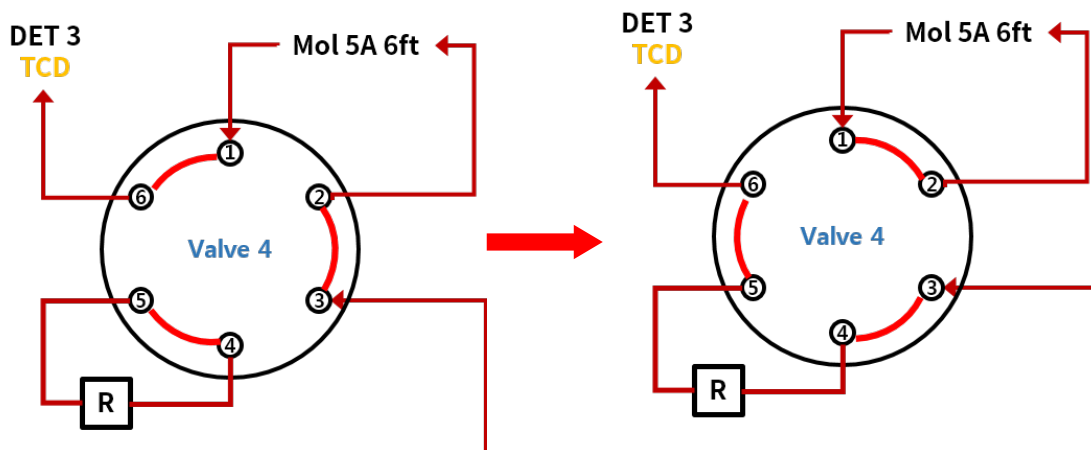
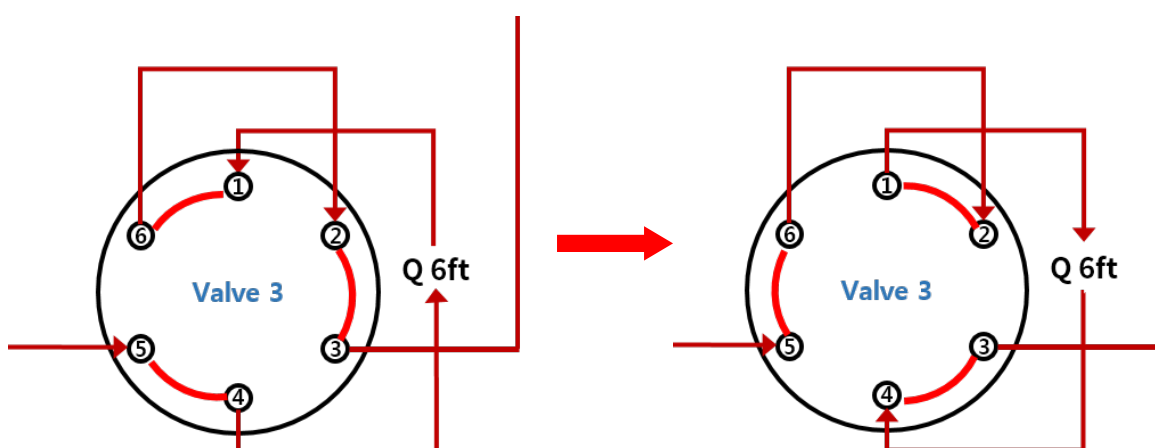
## Gas mixture analysis


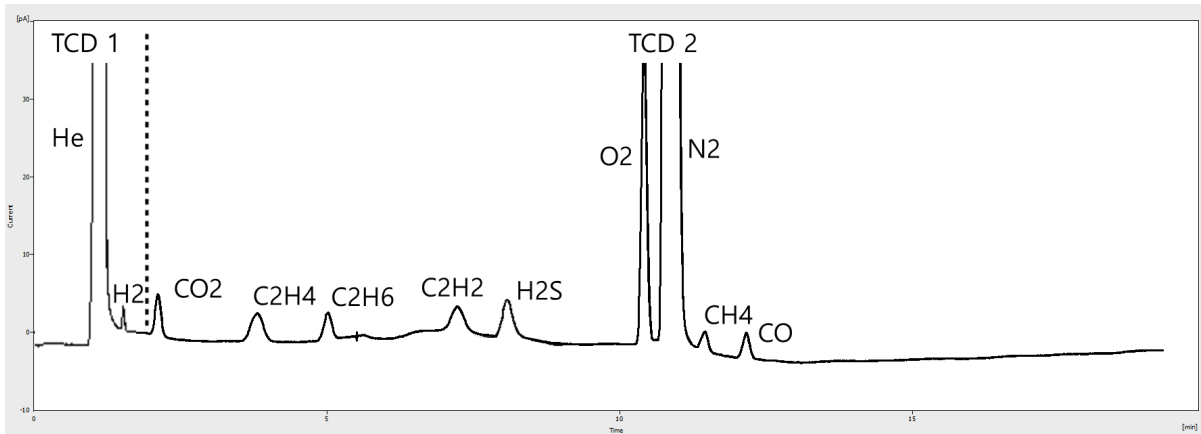
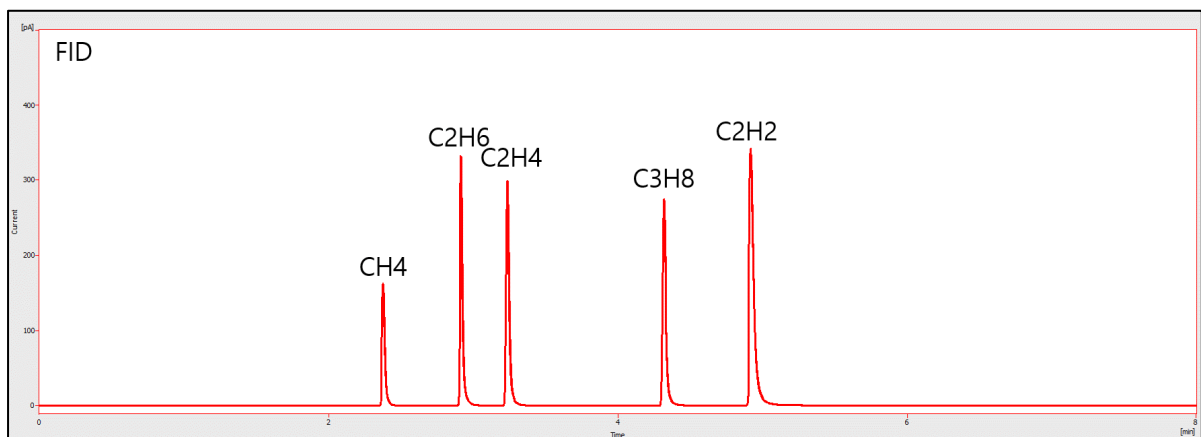
Fig 3. Chromatogram of Natural Gas Mixture

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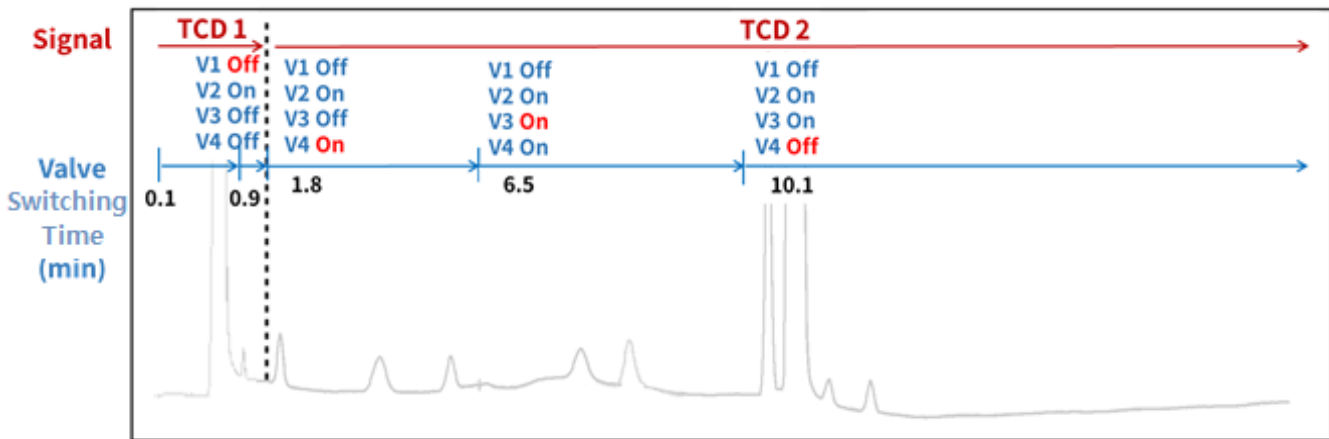## Conclusion

In this study, the determination of the chemical composition of natural gas mixture was conducted by ChroZen GC configuring 5 columns and 3 detectors with 4 valves according to ASTM D1945.

Also, it's easy to modify the valve configuration depending on the target gaseous sample and figure out the optimized valve switching time.

As the results, ChroZen GC verifies the right solution for superior resolution and sensitivity of components in natural gases within the range of composition shown in ASTM D1945.

## Reference

- ASTM D1945 Standard Test Method for Analysis of Natural Gas by Gas Chromatography

60, Anyangcheondong-ro, Dongan-gu,  
Anyang-si, Gyeonggi-do, 14042, Korea

TEL: +82-31-428-8700

FAX: +82-31-428-8787

E-mail: [export@youngincm.com](mailto:export@youngincm.com)

Homepage: [www.youngincm.com](http://www.youngincm.com)