


「应用说明」

甲苯、乙苯、二甲苯的ChroZen气相色谱分析

- GC 应用

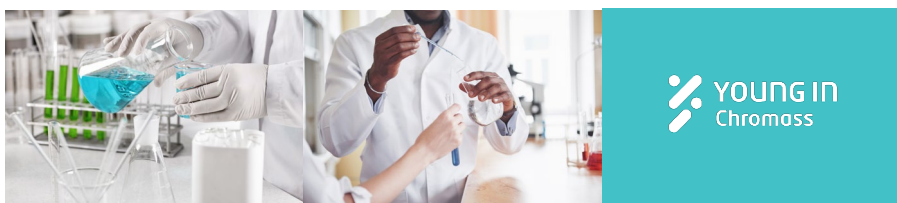

YOUNG IN
Chromass

分析条件

色谱柱: MEGA-5, 30 m x 0.32 mm x 0.25 μ m

柱温: 60°C(5 min), 40°C/min, 300°C(14 min)

载气: N₂


流速: 3 mL/min (20:1)

进样口: Capillary, 280°C

检测器: FID, 300°C

进样量: 1 μ L

结果

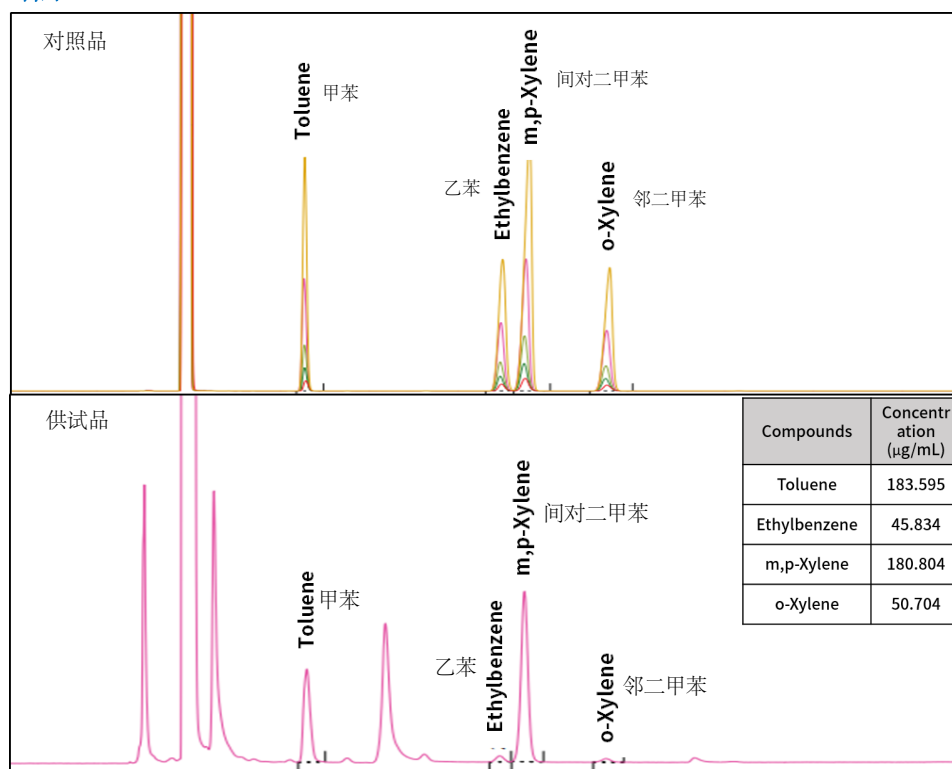

校准曲线

5 points – 25, 50, 100, 250, 500 µg/mL

