


初榨橄榄油中的甲酯、乙酯和蜡类

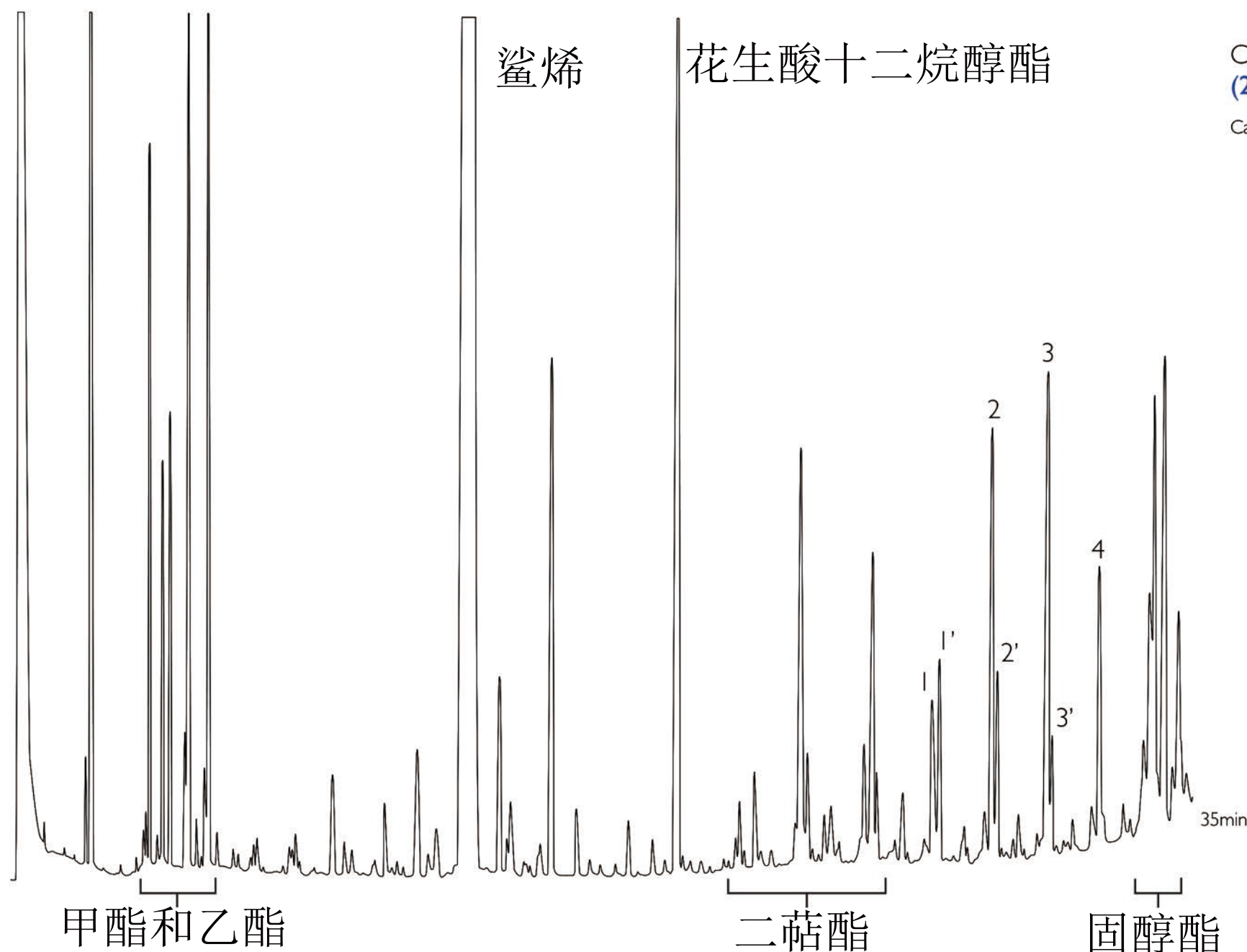

Column: **MEGA-SE52** - 0.32mm, 0.10 μ m, 10 + 2m
(2m of integrated built-in deactivated Retention Gap)

Catalog Code: GAP-SE52-032-010-10-2

蜡类分析的最优条件:

进样口: 冷柱上进样, 1 μ L正庚烷或者异辛烷溶液

检测器: FID 350 $^{\circ}$ C

柱温箱程序: 80 $^{\circ}$ C保持分钟后以20 $^{\circ}$ C

每分钟的升温速率升至220 $^{\circ}$ C, 然后以5 $^{\circ}$ C每分钟
升温至335 $^{\circ}$ C后恒温

载气: 氦气或者氢气, 选择最优线性的载气速度。

1+1':C40酯类

2+2':C42酯类

3+3':C44酯类

4+4':C46酯类

Acknowledgments: many thanks to **Prof. C. Mariani**, Stazione Sperimentale Oli e Grassi (SSOG), Milano, Italy.