


初榨橄榄油中的甲酯、乙酯和蜡类

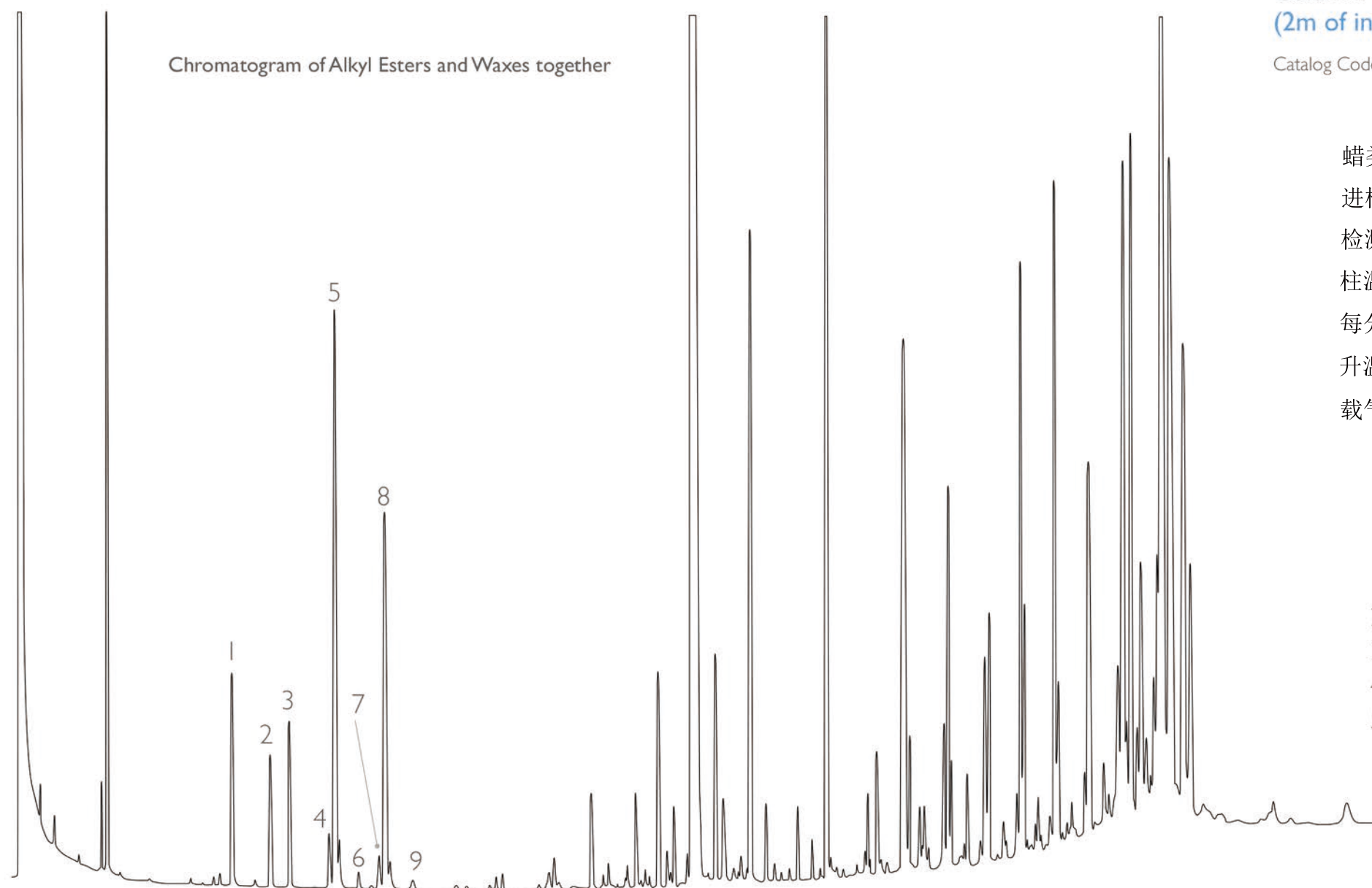

Column: MEGA-SE52 - 0.32mm, 0.10 μ m, 10 + 2m
(2m of integrated built-in deactivated Retention Gap)

Catalog Code: GAP-SE52-032-010-10-2

蜡类分析的最优条件:

进样口: 冷柱上进样, 1 μ L正庚烷或者异辛烷溶液

检测器: FID 350 $^{\circ}$ C

柱温箱程序: 80 $^{\circ}$ C保持分钟后以20 $^{\circ}$ C

每分钟的升温速率升至220 $^{\circ}$ C, 然后以5 $^{\circ}$ C每分钟
升温至335 $^{\circ}$ C后恒温

载气: 氮气或者氢气, 选择最优线性的载气速度。

- | | |
|------------------|-------------|
| 1. Me C16 | 6. Me C18 |
| 2. Et C16 | 7. Et C18:2 |
| 3. Me C17 (I.S.) | 8. Et C18:1 |
| 4. Me C18:2 | 9. Et C18 |
| 5. Me C18:1 | |

Acknowledgments: many thanks to Prof. C. Mariani, Stazione Sperimentale Oli e Grassi (SSOG), Milano, Italy.