

甘油三酯-榛子油

on MEGA-LAP Column

Column: **MEGA-LAP - 0.32mm, 0.10µm, 25m**

Catalog Code: C-LAP-032-010-25

MEGA-LAP (Lipid Analysis Phase) is a dedicated stationary phase optimized for Lipids, Sterols and saturated and unsaturated Triglycerides separations, stable over 370°C.


条件:

进样口: 分流 300°C, 1:80分流比, 1.0µL进样量

检测器: FID 300°C

柱温箱程序: 120-330°C, 5°C/min

载气: 氢气, 2mL/min

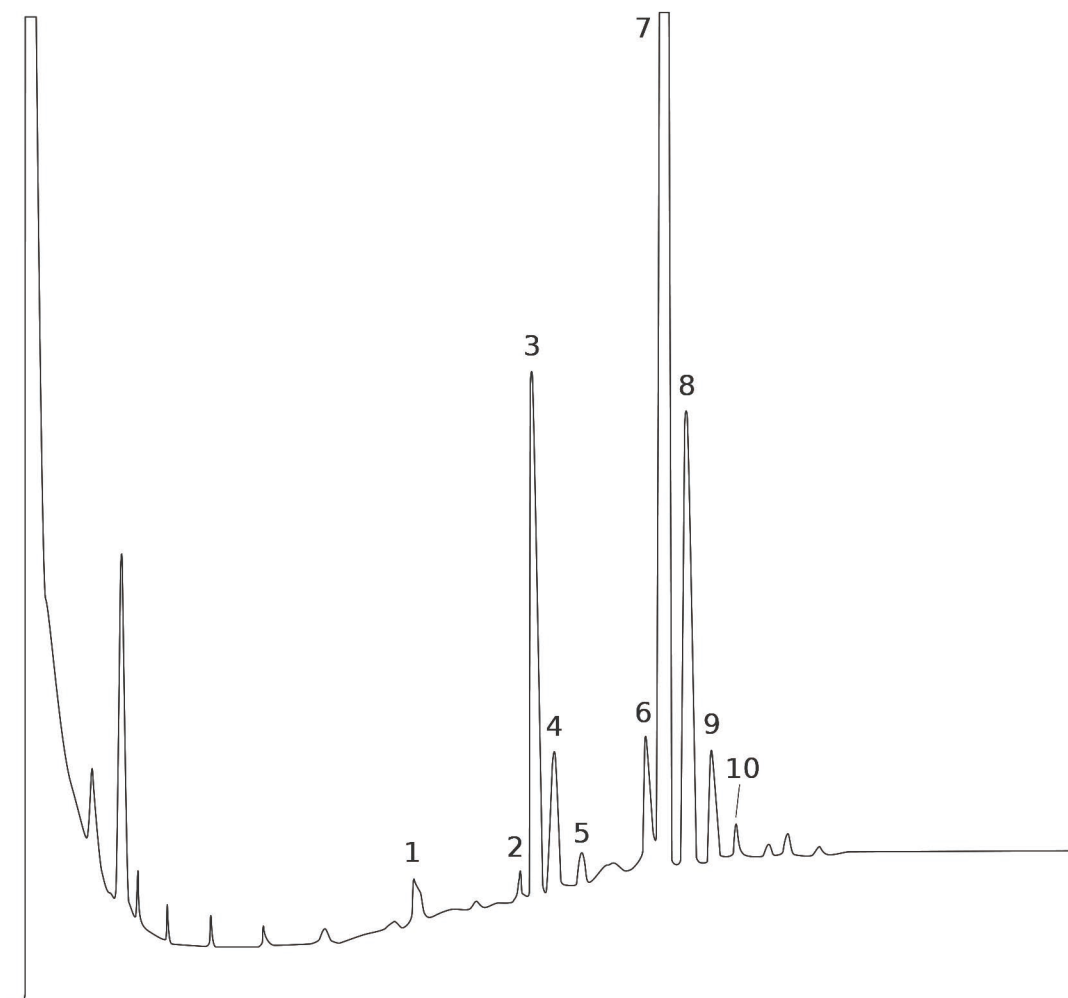

Peaks:

1. PPO
2. POS
3. POO
4. PLO + OOPo
5. PLL + PoOI
6. 8OO
7. OOO
8. OOL
9. OOL
10. LLL

Acknowledgement: Dott. C. Mariani - Stazione Sperimentale Oli e Grassi, Via Giuseppe Colombo 79 - Milano - Italy.