


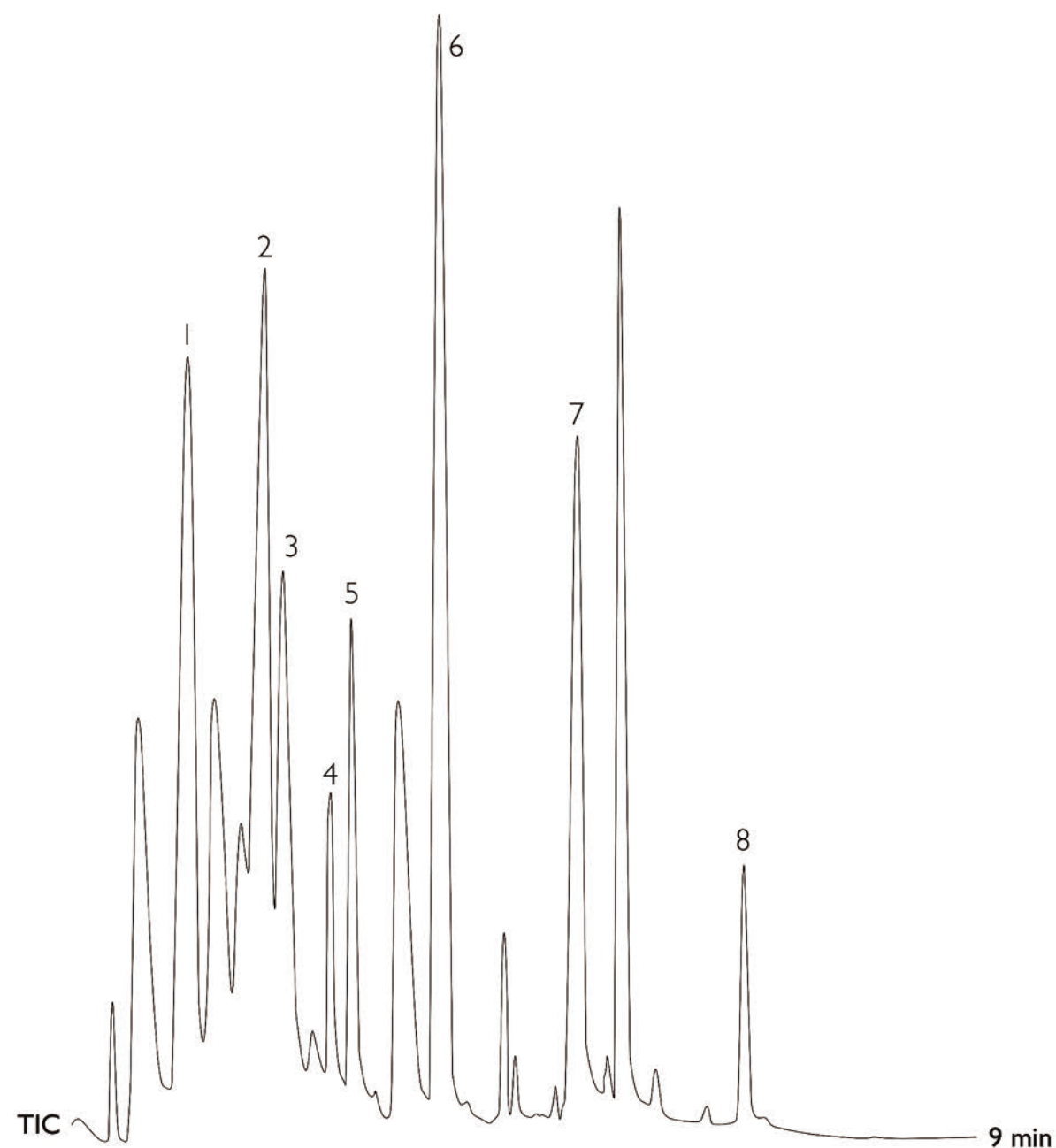

Analysis carried out with a DANI Master GC coupled with DANI Master TOF MS detector.

Column: **MEGA-5 FAST - 0.10mm, 0.10 μ m, 10m**
Catalog Code: F-5-010-010-10

条件:

进样口: 分流 280 $^{\circ}$ C, 1 μ L进样量

检测器: TOF-MS 系统

柱温箱程序: 50-300 $^{\circ}$ C (2min), 35 $^{\circ}$ C/min

载气: 氦气, 0.6mL/min

1	4.450 min	Phenacetin	非那西汀
2	4.900 min	Caffeine	咖啡因
3	5.000 min	Lidocaine	利多卡因
4	5.270 min	Levamisole	左旋咪唑
5	5.400 min	Procaine	普鲁卡因
6	5.920 min	Cocaine	可卡因
7	6.720 min	Ethylmorphine	乙基吗啡
8	7.690 min	Diltiazem	地尔硫卓