

- 1.丙三醇
- 2.丁三醇
- 3.脂肪酸(16) 甲酯
- 4.十六碳酸甲酯
- 5.单酰甘油
- 6.甘油单亚麻
- 7.单酰基甘油+亚油
- 8.十八烷酰
- 9.菜油甾醇TMS
- 10.豆甾醇TMS
- 11.β-谷甾醇
- 12.甘油三癸酸酯
- 13.甘油二油酸酯
- 14.β-谷甾醇葡萄糖苷TMS

Column: **MEGA-BIODIESEL 105** - 0.32mm, 0.10μm, 15m

Catalog Code: C-BIODIESEL105-032-010-15

条件:

进样口: PTV进样口, 温度100℃-370℃


检测器: FID 380℃

柱温箱程序: 50℃保持1min, 15℃/min, 升至180℃,
7℃/min, 升至230℃, 30℃/min, 升至370℃并保持15min

载气: 氮气, 3mL/min

Sample:

Biodiesel sample MSTFA derivatized, additioned with Butanetriol and Tricaprin (internal standards).

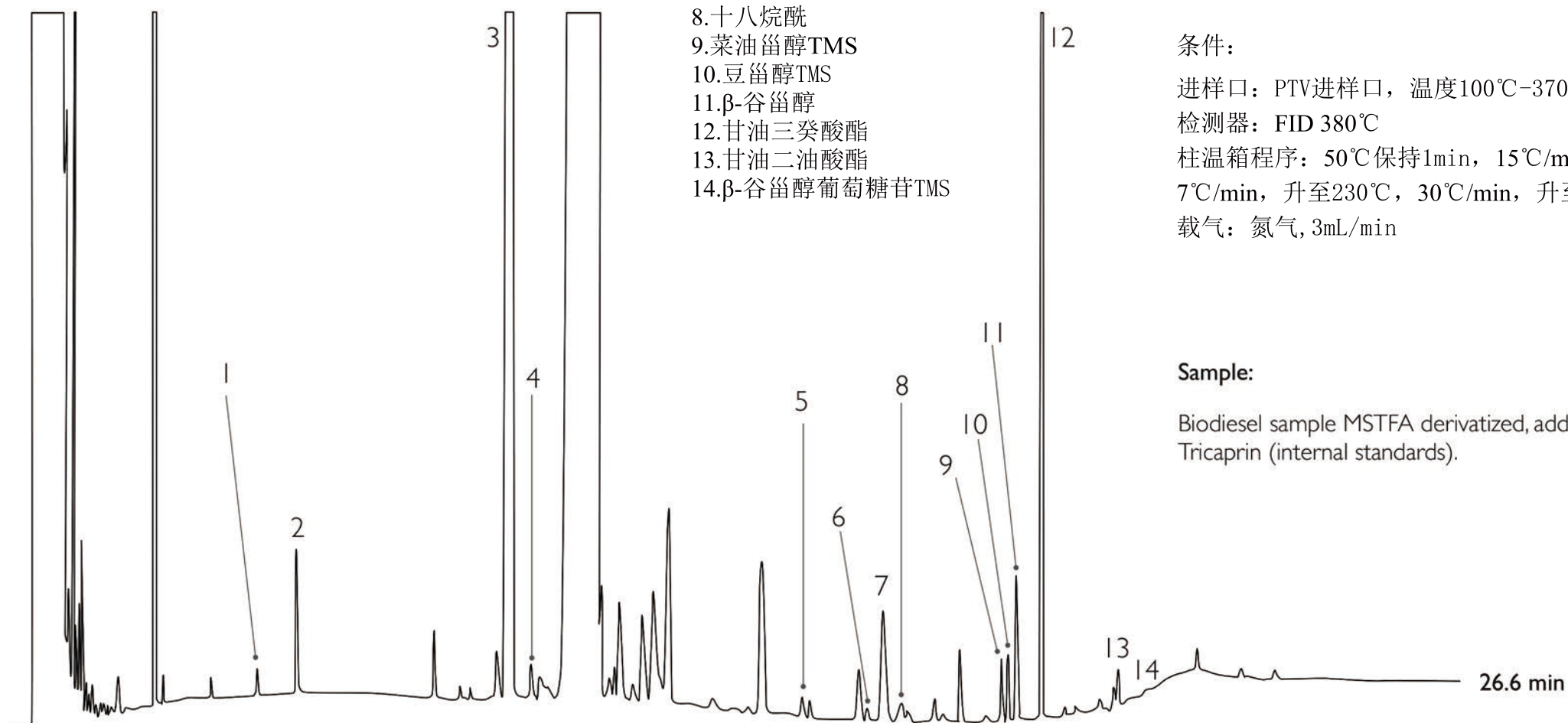

Acknowledgments: analysis carried out by **Lic. Rodolfo A. Cabrera** (Unitec Bio S.A., Santa Fe - Argentina). **Lab Solutions SRL** (Buenos Aires - Argentina - www.labsolutions.com.ar) is our official distributor and partner for Argentina.



Lab Solutions



MEGA®

www.mega.mi.it