

食品包装溶剂残留

on MEGA-VOC 1 & MEGA-VOC 2 dual system

MEGA-VOC 1 & 2 kit available for single injector port, dual detector GC system.

Contact us for more information.

条件:

进样口: 分流 250°C, 用微量进样针进样0.2μL,

150mL/min分流比

检测器: FID 250°C

柱温箱程序: 30-200°C, 35°C/min

载气: 氦气, 60kPa

峰:

- 1.乙醇
- 2.丙酮
- 3.2-丙醇
- 4.1-丙醇
- 5.丁酮
- 6.乙酸乙酯
- 7.环己烷
- 8.异丁醇
- 9.乙酸异丙酯
- 10.1-丁醇
- 11.甲氧基丙醇
- 12.乙酸丙酯
- 13.甲基异丁基酮
- 14.乙氧基丙醇
- 15.甲苯
- 16.乙酸丁酯

